Temperature

Controlled Shaking

Leading Safety Standards

Superior Ease of Use

Reduced Cost of Ownership

Operating Manual

Page 2

Test Tube Shaker Hei-MIX Multi Reax

Operating manual must be read before initial start-up.

Please follow the safety instructions provided.

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About this Document	4
Version and variants	4
About this manual	4
Symbols and keywords.....	5
Safety Instructions	6
General safety instructions.....	6
EU Declaration of Conformity.....	6
Intended use.....	6
Installation / electrical safety	7
Qualifications of employees.....	7
Obligations of the operator	7
Installation site	7
Modifications to the device.....	8
Employee safety	8
California Residents	8
Safety during use	9
Disposal	9
Device Description	10
Device overview	10
Overall view.....	10
Control panel	10
Start-Up	11
Setting up device	11
Set up.....	11
Power cord	11
Connect/disconnect power cord	11
Switch unit on and off	12
Carousel attachment.....	13
Attach vessels	13
Insert vessels.....	14
Operation	15
Operation and controls	15
Continuous shaking.....	16
Timer controlled shaking.....	16
Repetition	16
Edit time during shaking operation	16

Cleaning and Maintenance	17
Cleaning.....	17
Maintenance	17
Troubleshooting	18
Possible errors and how to resolve	18
Assembly	19
Electrical connection	19
Exchange carousel attachment	20
Disassembly and Storage	21
Dismantling, transportation and storage	21
Dismantling	21
Transportation and storage	21
Accessories and Spare Parts	22
Scope of delivery	22
Accessories.....	22
Attachments	23
Technical data	23
Service	26
Contact / Technical Service	26
Warranty	27
Confirmation of condition.....	28

Copyright

This operating manual is protected by copyright and is exclusively intended for the internal use of the purchaser.

Handing over this operating manual to third parties, reproducing it in any type or form – in part or in full – and utilizing and/or communicating its contents is prohibited without prior written approval from the manufacturer.

Any violations will result in an obligation to provide compensation for damages. Other claims remain unaffected.

➤ Version and variants

Version

This manual describes the function, operation and maintenance of the Hei-MIX Multi Reax test tube shaker.

Version	Alteration date
1.2	12/2019

Variants

The device is available in different variants. Certain features or functions are only available in specific product variants. These variants are described in this operating manual.

➤ About this manual

This operating manual is an integral part of the device described here.

- Carefully read through this operating manual and observe all of the safety instructions and warnings notices.
- Ensure that every user has carefully read the operating manual before using the device for the first time.
- Store this operating manual so that it is accessible to all users at all times.
- Pass the operating manual on to any subsequent owner.

The current version of this operating manual can also be found at:
[https://heidolph-instruments.com/en/service/downloads/operation-manuals#Shakers & Mixers \(Multi Reax\)](https://heidolph-instruments.com/en/service/downloads/operation-manuals#Shakers & Mixers (Multi Reax))

In addition, also observe the local regulations at the installation site.

➤ Symbols and keywords

Standardized symbols and signal words are used in this operating manual to warn against any dangers and provide important instructions. These instructions must be strictly observed to avoid accidents and damage.

The following terms and basic symbols are used:

Symbol	Additional signal words / explanation
Warning signs 	<p>The yellow triangle indicates hazardous situations. It is used in combination with the following signal words:</p> <p>DANGER: Indicates a hazardous situation which, if not avoided, will result in serious injury or death.</p> <p>WARNING: Indicates a hazardous situation which, if not avoided, may result in serious injury or death.</p> <p>CAUTION: Indicates a hazardous situation which, if not avoided, may result in property damage and minor or moderate injury.</p>
Prohibitory signs 	<p>Prohibited:</p> <p>The red circle indicates a situation that should be avoided under all circumstances and which, if not avoided, may result in serious injury or death.</p>
Mandatory signs 	<p>Must be observed:</p> <p>The blue circle indicates important information. Please observe this information to avoid any property damage.</p>

Other symbols used:

Symbol	Description
➔	Handling instruction, action required
✓	Result of action
▪	List of information
▪	
▪	
a.	List of variants
b.	
c.	

➤ General safety instructions

EU Declaration of Conformity

This device complies with the following EC-Directives*:

- 2006/42/EC Machinery Directive
- 2014/30/EU Electromagnetic Compatibility Directive

*Also see the attachment "EU Declaration of Conformity".

The device has been manufactured according to state-of-the-art technology and in compliance with recognized safety regulations. However, risks may still arise during installation, operation and maintenance.

→ Please ensure that the operating manual is available at all times.

The device may only be used under the following conditions:

- Only operate the device if it is in a technically perfect condition.
- Only operate the device if you are properly aware of the risks and required safety measures.
- Only operate the device in accordance with the instructions given in this operating manual.
- If any information is ambiguous or missing, ask your superior or contact the manufacturer.
- Do not operate anything on the device without authorization.
- Only use the device in accordance with its intended use.

Intended use

The device may only be operated by authorized personnel.

The device is suitable for the following uses:

- Shaking
- Mixing
- Emulsifying
- Suspending
- Dissolving

The device may be operated in research laboratories, other laboratories and production facilities in the following sectors:

- Chemistry
- Pharmacy
- Biology
- Environmental analytics
- Basic research
- Similar research laboratories

Unintended use/Reasonable foreseeable misuse

Any use which deviates from the device's intended use is considered to be improper. The manufacturer is not liable for damage that occurs as a result. The risk is borne by the operator alone.

When used in medical and food sectors the operator alone is responsible for compliance with the applicable regulations and standards.

Installation / electrical safety

- The device may only be connected up if the available power supply voltage corresponds to the information stated on the rating plate for the device.
- The power supply connection must be easily accessible at all times.
- An RCD circuit breaker (residual current circuit breaker) must be fitted to the electrical system in the building so that the device is disconnected from the power circuit in the event of a fault.
- Repairs may only be performed by electricians authorised by Heidolph Instruments.
- Never operate the device with a damaged power cord.
- Always turn the device OFF and disconnect the power cord before carrying out any maintenance or repair work.

Qualifications of employees

- The device may only be operated by trained personnel.
- The device may only be operated by persons who have been instructed and supervised in its proper use by trained specialist personnel.
- The device may only be operated by specialist personnel who are above the legal minimum age.
- Other persons may only work on the device under the constant supervision of experienced and trained specialist personnel.
- This operating manual must be read and understood by all persons working with the device.
- Personnel must receive safety training that ensures responsible and safe working practices.

➤ Obligations of the operator

Installation site

- The device must be installed in a suitable location.
- The device must be installed on a firm and stable surface.
- Ensure that the device and all of its components are easily accessible at all times.
- Maintain a sufficient safety clearance to, in particular, any moving and/or hot equipment components.
- It is not permitted to place or store any objects such as accessories, tools or chemicals within this clearance area during operation.
- All screw connections must be securely tightened.
- It is not permitted to operate the device near to highly flammable or explosive substances.
- Operate the device in conjunction with an extractor hood, if working with potential harmful media (see DIN EN 14175 and DIN 12924).

Safety Instructions

- The device is designed for indoor use in dry spaces only and under the following ambient conditions:

Ambient temperature	5 - 31 °C at 80 % relative humidity
	32 - 40 °C decreasing linearly to a maximum 50 % relative humidity
Installation altitude	0 - 2,000 m above sea level
Contamination level	2
Overvoltage category	II
Permissible supply deviations	± 10 %

- If the device is operated in corrosive atmospheres, the service life of the device will decrease based on the concentration, duration and frequency of the exposure to the corrosive atmosphere e.g. concentrated hydrochloric acid (HCl).

Modifications to the device

- It is not permitted to make any unauthorised modifications or changes to the device.
- Do not attach or install any parts that have not been approved by the manufacturer.
- Unauthorised modifications or changes will void the EC Declaration of Conformity for the device and operation of the device will no longer be permitted.
- The manufacturer is not liable for any damage, dangers or injuries that result from unauthorised modifications and changes or due to the non-observance of the instructions in this manual.

Employee safety

- Ensure that the device is only operated by qualified specialist personnel and trained employees.
- Observe the following instructions to avoid any personal injuries and property damage:
 - Laboratory regulations
 - Accident prevention regulations
 - Hazardous Substance Act
 - Other generally accepted occupational health and safety regulations
 - Local regulations

California Residents

- Important information for California residents regarding Prop 65. Please visit www.P65Warnings.ca.gov for more information.

➤ Safety during use

- Wear the appropriate protective clothing when working on the device (protective glasses and, if necessary, safety gloves).
- Do not use the device in potentially explosive areas. The device is not protected against explosion. There is no explosion or ATEX protection available.
- Do not carry out work with naked flames in the vicinity of the device (risk of explosion).
- Do not operate any devices in the vicinity which may exhibit electromagnetic fields in the frequency range 9×10^3 Hz to 3×10^{11} Hz.
- Do not operate or assemble devices in the vicinity which are emission or radiation sources (electromagnetic radiation) for the frequency range 3×10^{11} Hz to 3×10^{15} Hz (in the optical spectral range wavelengths from $1,000\mu$ to 0.1μ).
- Do not operate or assemble appliances in the vicinity of the device which constitute emission or radiation sources for ionizing radiation or in the ultrasonic range.
- Do not operate the device where adiabatic compression or shock waves might occur (shock wave combustion).
- Do not use substances, where the energy input due to mixing might pose a hazard.
- Avoid putting pressure on the display when you are not operating the device.
- Do not spill liquids over the device or any parts of it.
- Remove any accidentally spilled liquids immediately.
- Eliminate errors immediately.
- Do not use abrasive material to clean the surface. Only wipe with damp cloth.
- Always switch the device OFF after use.

➤ Disposal

- Check the device components for hazardous substances and solvents.
- Clean all components before disposal.
- Dispose of the device in accordance with the relevant national regulations.
- Dispose of the packaging material in accordance with the appropriate national regulations.


➤ Device overview

Overall view

Control panel

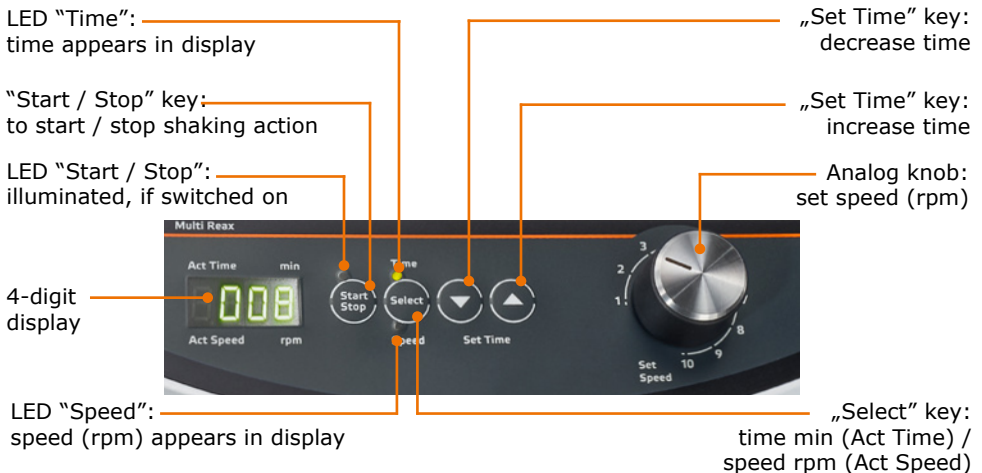➤ Setting up device

Warning:
Risk of slipping device

As a result of vibrations generated during operation the device might slide off the table top.

The surface must be smooth and clean and should be properly leveled.

Set up

- Locate the shaker on a stable, horizontal surface.
- Clean surface and feet with a damp cloth and ethanol regularly.

➤ Power cord

Connect/disconnect power cord

Connect plug to device

The power cord comes with a three-wire plug and a recess on the bottom side.

Unit plug

Power cord

The socket is located on the back side of the unit.

- ✓ The device is switched OFF.
- Connect plug to the socket with the recess facing down.
- Push plug up against socket until securely attached.
- Connect plug to power socket.

Unit socket

Disconnect power cord

- Disconnect plug from power socket first.
- Then disconnect plug from socket on unit.

➤ Switch unit on and off

The ON/OFF switch is situated on the right side of the device beneath the control panel.

Switch unit on

- Press the ON/OFF switch on the upper side.
- ✓ The unit is switched on.

Switch unit off

- Press the ON/OFF switch on the lower side.
- ✓ The unit powers off.

Switch OFF

Switch ON

➤ Carousel attachment

Attach vessels

Caution:
Risk of breakage and spilling

If the rubber tension ring is damaged your vessels may fall out of the attachment and spill.

Control the rubber tension ring regularly. Replace it eventually.

Vessel sizes and load

- Carousel attachment for 26 test tubes
- round vessels, Ø 10 – 16 mm
- max. length 160 mm
- Carousel attachment for 12 test tubes
- round vessels, Ø 16 – 32 mm
- max. length 120 mm.
- Maximum load for device
- 1.5 kg

Attach vessels

- ➔ Plug vessels or tubes into holding pockets against preload of a rubber tension ring.
- ✓ This rubber tension ring is going to safely secure the tubes in the holding position.

Carousel attachment

Rubber tension ring

Holding pockets

- ➔ To ease installation, lift rubber tension ring by hand before introducing vessels.


Insert vessels

Caution:
Risk of breakage and spillage

If vessels are not inserted properly they may fall out of the attachment and spill.


Refer to recommendations below when inserting vessels.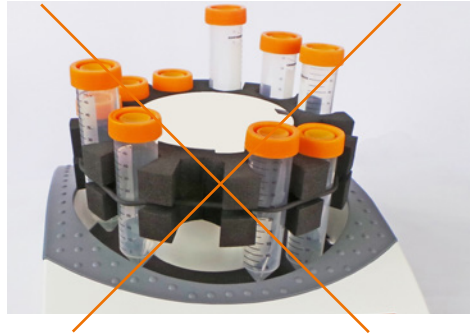


- Ensure that rubber tension ring holds the center of the vessel.

- Distribute vessels around evenly.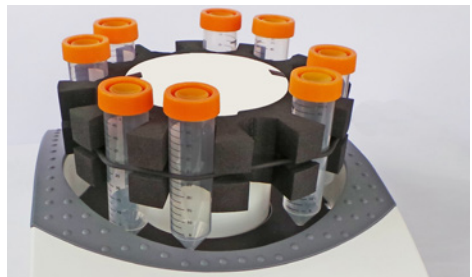

- **Never** attach different vessel sizes randomly. The smaller diameter vessel will not be securely fastened!
- Form groups of vessels with same diameter and leave at least one space between groups open.

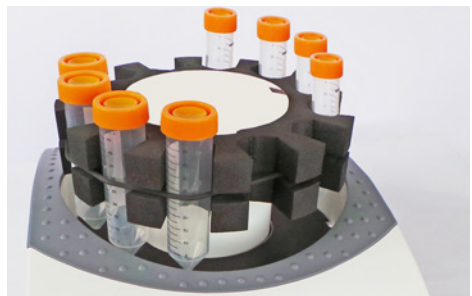➤ Operation and controls

Caution:
Risk of breakage and spillage

If the vessels are not attached securely they may fall out and spill the sample.

Prior to switching the unit on ensure that all vessels are securely fastened.

Warning:
Risk of poisoning

Open vessels and too high vibration frequencies may result in samples splashing.

Close all vessels tight especially if sample is dangerous or toxic.

Adjust vibration frequency step by step until you have reached the required rpm settings..

Wear safety glasses and especially with dangerous and toxic samples appropriate protective clothing.

Warning:
Risk of injury!

High speeds in combination with high loads and/or a tall assembly may lead to an unstable device. As result the device may start moving and slide off the table top.

When working with with high loads and/or a tall assembly, increase the speed with care and watch for the devices stability. Ensure the device is installed in such a way as it can not collide with the surrounding environment or slip off its work surface.

Before switching the unit on ensure that no vessel can strike against the housing.

Depending on the quantity and the load of vessels, certain speed levels may create a resonance. Increase or decrease speed by a few rpm until the resonance disappears.

A power cut will stop the unit.
With power return the device is switched on automatically. Shaking must be started manually.

Continuous shaking

- Turn device ON with ON/OFF switch.
- Set time with ↑ and ↓ keys to 000 (depress ↑ and ↓ together).
- Start continuous operation by hitting the "Start / Stop" key.
- ✓ LED "Start" and "Speed" turn on, shaking frequency is displayed (4-digit display).
- Set shaking frequency with knob.
- Stop shaking Hitting "Start / Stop" key will stop.

Timer controlled shaking

- Set time between 0 and 999 minutes with Set Time ↑ or Set Time ↓ keys.
- ✓ Short hits on the keys change time setting slowly, whereas depressing keys is going to change time rapidly.
- ✓ Depressing the Set Time ↑ and Set Time ↓ keys resets the timer to 000.
- Start shaking by pressing the "Start / Stop" key at the frequency selected with knob and for the time interval set before.
- ✓ LED "Start" turns ON.
- ✓ Time lapse appears in a 4-digit display. LED "Time" turns ON.
- Discontinue current shaking action by pressing the "Start / Stop" key.
- ✓ Actual time won't change.
- Continue shaking by hitting the "Start / Stop" key.
- ✓ In case of a power supply failure or interruption (item turned OFF), the time set is lost.
- Change display from time to shaking frequency and vice versa by hitting the "Select" key during timer-controlled operation.
- ✓ LED "Speed" turns ON. Timer LED flashes to indicate timer operation mode.
- ✓ LED "Time" also flashes, when timer-controlled operation was discontinued with the "Start / Stop" key.
- ✓ As soon as time has elapsed shaking stops and an acoustic signal will be audible.

Repetition

Repeating last timer controlled operation does not require new time setting, as long as you did not touch the master switch.

Edit time during shaking operation

- Time setting may be changed during shaking, while timer is running and even when timer controlled operation was discontinued.
- The timer will keep the initial value in its memory for an eventual repetition function.

> Cleaning

All surfaces of the unit allow for cleaning with a damp cloth and if required a mild soap lotion.

Warning:
Risk of electric shock!

If the device is still connected to power during cleaning, you might get in contact with live parts due to accidentally entering liquids.

Prior to cleaning:

- ensure that the device is switched OFF
 - ensure that the plug is pulled from the electric socket.
-

Caution:
Damage to the device

Accidentally entering liquids will cause the device to fail .

Only wipe with a damp cloth. Never use a soaked cloth.

Attention:
Risk of damage to the surfaces

All of the surfaces may be damaged by improper cleaning.

Never use:

- Chlorine bleach or any cleaning agent containing chlorine
 - Solvent-bearing agents (e.g. acetone)
 - Ammonia
 - Abrasive cleaning agents such as scouring agents, cleaning wool or agents with metallic particles
-

As a result of vibrations generated during operation the device might slide off the table top.

To ensure a stable position clean surface and feet with a damp cloth and ethanol regularly.

> Maintenance

The unit is maintenance-free. Any necessary repair must be performed by an authorized Heidolph distributor.

Please contact Heidolph Instruments or your local Heidolph Instruments distributor.

Troubleshooting

➤ Possible errors and how to resolve

Malfunctions	Possible reason	Troubleshooting
Unit won't shake	No power	<ul style="list-style-type: none"> Check size and shape of plug and compatibility to your electrical socket. Check circuit breakers
Unit stops shaking	Pre-set timer elapsed	Restart timer, if necessary
	Thermal motor circuit breaker triggered by motor overload.	Wait 20 minutes, decrease overall load
	<ul style="list-style-type: none"> Mechanical parts broken Electronic failure 	Contact your local Heidolph Instruments distributor

If you experience a malfunction which can not be resolved, please contact your authorized Heidolph Instruments distributor immediately.

➤ Electrical connection

Repairs to the device must **only be carried out by a qualified electrician approved by Heidolph Instruments**. Improper repairs could result in serious dangers. Please contact your specialist dealer if the device needs to be repaired.

Warning: **Risk of electric shock**

The device must be connected to an earthed power socket.

The rated voltage for the device and the supply voltage must be the same. The rated voltage for the device can be found on the rating plate on the rear of the device.

The device is earthed via the supplied power cord when delivered. If the original plug is replaced, it is essential that the new plug is connected to a protective earth!

Rating plate

Colour coding for the electrical connection:

Europe

**GREEN/
YELLOW** PE: Protective earth
(**earth**)

BLUE N: Neutral conductor

BROWN P: Phase

North America

GREEN PE: Protective earth
(**earth**)

WHITE N: Neutral conductor

BLACK P: Phase

The device is connected to the power supply with the supplied power cord. The device socket is located on the rear of the device.

If you want to use the device in a country with a different type of plug:

- The supplied plug may only be replaced by a qualified electrician.
- If you want to use an adapter, it must be earthed and approved in accordance with local regulations.
- The cable, adapter and plug must correspond at least to the power stated on the rating plate.


➤ Exchange carousel attachment

The item arrives with the carousel attachment for 26 test tubes installed. You may replace it with the carousel attachment for 12 test tubes at ease.

- Remove all vessels from carousel attachment.


- Remove carousel attachment by pulling upwards cautiously off the two fixing noses with both hands.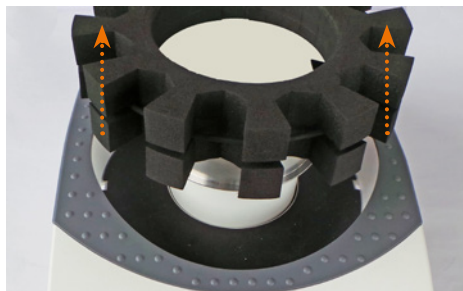


Be sure not to damage the attachment by the fixing noses. The fixing noses secure the carousel attachment on the shaker.

- Carefully slip carousel attachment over the two fixing noses in place.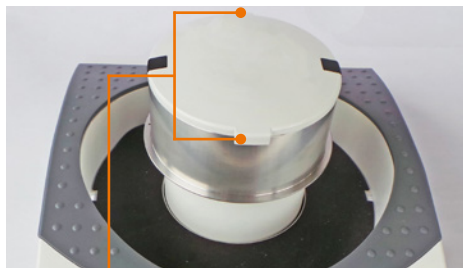

Haltenasen

- Check rubber tension ring for proper seat. This ring secures your vessels in place.
- Seat ring in slot of carousel if applicable.

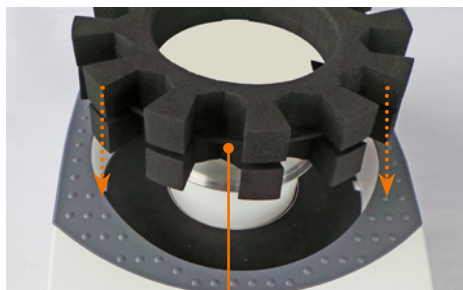Rubber tension ring

➤ Dismantling, transportation and storage

Dismantling

Warning:
Risk of injury!

While the device is plugged in you might be at harm if the device is accidentally switched to ON mode.

Prior to dismantling device ensure that the device is switched off and the plug is pulled from the electric socket.

- ✓ The device is switched off.
- Disconnect plug from power socket.
- Disconnect plug from socket on the device.
- Remove all vessels.
- Disassemble attachments.

Transportation and storage

- Store and transport the device and its components only if they were emptied, cleaned and dismantled.
- To carry the device wear safety gloves and lift the unit from below.
- Store and transport the device and parts of device in the original packing material or alternatively in an adequate container to prevent damage.
- Seal the package carefully against unauthorized or accidental opening.
- Store the device in a dry and frost-free place.

Improper storage and transportation may result in damages to the system and the mechanical components.

Avoid any kind of shocks during transportation.

➤ Scope of delivery

Component	Quantity	Product number
Hei-MIX Multi Reax*	1	545-10000-00
Carousel attachment for 26 test tubes, 10 - 16mm Ø	1	549-01600-00
Carousel attachment for 12 test tubes, 16 - 32mm Ø	1	549-03000-00
Operating Manual English/German	1	01-005-004-28
Warranty registration / Confirmation of condition	1	01-006-002-78
EC Declaration of Conformity	1	01-001-025-06
Power cord	1	depending on country

* The product number is valid for 230 V devices in the european community. Please contact Heidolph Instruments for further product numbers on country specific varieties.

➤ Accessories

Component	Quantity	Product number
Rubber tension ring for carousel attachment 26 tubes	1	23-08-06-01-75
Rubber tension ring for carousel attachment 12 tubes	1	23-08-06-01-76

Additional accessories are shown in the general catalog or on our website at www.heidolph-instruments.com

➤ Technical data

Multi Reax

Standard supply voltage	230 V/50 Hz or 115 V/60 Hz
Input power (W)	50
Protective class (IEC 61140)	1 ⚡
Protection class (IEC 60529)	IP 30
Sound pressure level (dB(A)) (based on IEC 61010)	< 45
Drive	asynchronous motor
Overheat protection	self-resetting
Rotation speed (rpm)	150 - 2,000
Rotation speed setting	digital
Motion	circular vibrating
Orbit (mm)	3
Load capacity (kg)	1.5
Operating mode	automatic or continuous
Weight (kg)	9.8
Dimensions (W x h x d)	270 x 172 x 410

China RoHS DECLARATION OF CONFORMITY

Heidolph Instruments GmbH & Co.KG has made reasonable efforts to ensure that hazardous materials and substances may not be used in its products.

In order to determine the concentration of hazardous substances in all homogeneous materials of the subassemblies, a "Product Conformity Assessment" (PCA) procedure was performed. As defined in GB/T 26572 the "Maximum Concentration Value" limits (MCV) apply to these restricted substances:

- Lead (Pb): 0.1%
- Mercury (Hg): 0.1%
- Cadmium (Cd): 0.01%
- Hexavalent chromium (Cr(VI)): 0.1%
- Polybrominated biphenyls (PBB): 0.1%
- Polybrominated diphenyl ether (PBDE): 0.1%

Environmental Friendly Use Period (EFUP)

EFUP defines the period in years during which the hazardous substances contained in electrical and electronic products will not leak or mutate under normal operating conditions. During normal use by the user such electrical and electronic products will not result in serious environmental pollution, cause serious bodily injury or damage to the user's assets.

The Environmental Friendly Use Period for Heidolph Instruments GmbH & Co.KG products is 25 years.

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This table is created according to SJ/T 11364-2014

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Note: Table applies to all products. Some of the components or parts listed above may not be part of the enclosed product.

- O: 表示该有毒有害物质在该部件所有均质材料中的含量均在GB/T 26572规定的限量要求以下。
O: Indicates that the above mentioned hazardous substance contained in all homogeneous materials of the part is below the required limit as defined in GB/T 26572.
- X: 表示该有毒有害物质至少在该部件某一均质材料中的含量超出GB/T 26572规定的限量要求。
X: Indicates that the above mentioned hazardous substance contained in at least one of the homogeneous materials of this part is above the required limit as defined in GB/T 26572.

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Apart from the disclosures in the above table, the subassemblies are not intentionally manufactured or formulated with lead (Pb), mercury (Hg), cadmium (Cd), hexavalent chromium (Cr(VI)), polybrominated biphenyls (PBB), and polybrominated diphenyl ethers (PBDE).

Products manufactured by Heidolph Instruments GmbH & Co.KG may enter into further devices or can be used together with other appliances.

With these products and appliances in particular, Heidolph Instruments GmbH & Co.KG will not take responsibility for the EFUP of those products and appliances.

Place, date Schwabach, 19.09. 2019

Wolfgang Jaenicke
Chief Executive Officer CEO

Marcell Sarré
Vice President
Quality Management &
Technical Service

➤ Contact / Technical Service

Questions / Repair work

If any aspect of installation, operation or maintenance remains unanswered in the present manual, please get in touch with the following address.

For repairs please call Heidolph Instruments or your local authorized Heidolph Instruments distributor.

Warning! Danger of poisoning!

Contaminated devices can lead to severe injury or death of our employees!

When shipping items for repair that may have been contaminated by hazardous substances, please:

- advise exact substance
- take adequate protective action towards our parts receiving and service personnel
- mark the pack in accordance with Ordinance on Hazardous Substances

A "Confirmation of Condition" form can be found at the end of this operating manual.

→ Prior to shipping the device for repair, complete a copy of this form and submit it in advance.

Contact details:

Heidolph Instruments Germany

Heidolph Instruments GmbH & Co. KG
Technical Service
Walpersdorfer Str. 12
D-91126 Schwabach / Germany

Tel.: +49 - 9122 - 9920-74
Fax: +49 - 9122 - 9920-84

E-Mail: service@heidolph.de
www.heidolph.com

Heidolph Instruments North America

Phone: 1-866-650-9604
E-mail: service@heidolph.com
www.heidolphNA.com

Heidolph Instruments United Kingdom

Phone: 01799 - 5133-20
E-mail: service@radleys.co.uk
www.heidolph-instruments.co.uk

All other countries

You will find contact details of your local Heidolph distributor at
www.heidolph.com

➤ Warranty

Heidolph Instruments provides a 3 year warranty for the products described here (excluding glass and wearing parts) if you register using the warranty card enclosed or online (www.heidolph.com). The warranty period begins with the date of purchase. The serial number is also valid without registration.

This warranty covers material and manufacturing defects.

In the event of a material or manufacturing defect, the device shall be repaired or replaced free of charge under the terms of the warranty.

Heidolph Instruments shall not assume liability for any damage incurred as a result of improper handling or transport.

Warranty claim?

- ➔ Please notify Heidolph Instruments or your local Heidolph distributor if you wish to make a warranty claim.

➤ Confirmation of condition

→ In the case of repair, copy and complete the Confirmation of condition form and send it to Heidolph Instruments.

1. Details about the device

Product number

Serial number

Reason for repair

2. Has the device been cleaned, decontaminated/sterilized?

Yes

No

3. Is the device in a condition which does not represent any health threats for the staff of our service department?

Yes

No

If not, which substances has the device
come into contact with?

4. Legally binding declaration

The customer is aware of being legally liable to Heidolph Instruments
for any damages arising from incomplete and incorrect information.

Date

Signature

Company stamp

Please note

The shipper is responsible for the return of the goods in well packed
condition, suitable for the mode of transport.

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This is a translation of the original operating manual. We reserve the right to make any technical modifications without prior notification. This documentation is not subject to revision; the latest version of this documentation can be found on our homepage. The copyright for all texts and images is held by Heidolph Instruments GmbH & Co. KG.
© Heidolph Instruments GmbH & Co. KG