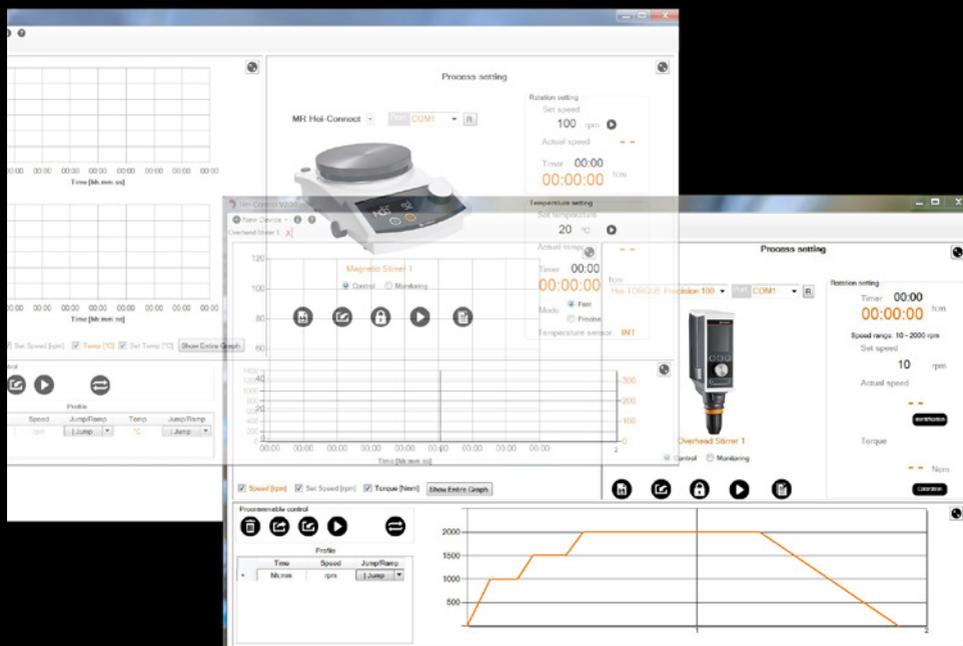
# PC controlled Stirring

Leading Safety Standards

Superior Ease of Use

Reduced Cost of Ownership

## Operating Manual

Page 2

Software Hei-CONTROL for overhead stirrer Hei-TORQUE and magnetic stirrer Hei-Connect

Operating manual must be read before initial start-up.

Please follow the safety instructions provided.

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<b>About this Document</b>	<b>4</b>
Version and variants .....	4
System requirements .....	4
About this manual .....	4
Terms and signs used.....	5
<b>Safety Instructions</b>	<b>6</b>
General safety instructions.....	6
Intended use.....	6
Operating company's obligations.....	6
Changes to the software.....	6
Personnel safety .....	6
Safety during use .....	7
Disposal .....	7
<b>Installation</b>	<b>8</b>
Installation .....	8
USB interface .....	8
Software Hei-Control.....	8
Manual Hei-Control .....	9
Installation adjustments .....	10
Updates .....	10
<b>Start-Up</b>	<b>11</b>
Interface .....	11
Start and end program .....	12
Link device and PC.....	12
Cut link between device and PC .....	13
<b>Overhead Stirrer Operation</b>	<b>14</b>
Process parameters .....	14
Menu structure.....	15
Register and remove several Overhead Stirrers.....	17
Operate or monitor stirrer.....	17
Symbols used.....	18
Control the Overhead Stirrer .....	19
Set stirring parameters.....	19
Start and Stop Stirring .....	20

Profile .....	21
Edit profile .....	21
Save profile, recall and reset .....	21
Delete profile .....	22
Start profile, interrupt, repeat, stop .....	22
Calibrate torque .....	23
Read out Data .....	23
Save and restore process data .....	23
Lock and unlock software .....	24

## **Magnetic Stirrer Operation** **25**

Process parameters .....	25
Menu structure.....	26
Register and remove several Magnetic Stirrers.....	28
Operate or monitor stirrers .....	28
Symbols used.....	29
Control the Magnetic Stirrer .....	30
Set stirring parameters.....	30
Start and stop stirring .....	31
Set heating parameters .....	32
Start and stop heating.....	33
Profile .....	34
Edit profile.....	34
Save profile, recall and reset .....	34
Delete profile .....	35
Start profile, interrupt, repeat, stop .....	35
Read out Data .....	36
Save and restore process data .....	37
Lock and unlock software .....	37

## **Troubleshooting** **38**

Malfunctions and clear up .....	38
---------------------------------	----

## **Service** **41**

Contact / Technical Service .....	41
-----------------------------------	----

### > Version and variants

#### Version

This manual describes the function and operation of the software program Hei-CONTROL for Hei-TORQUE Precision Overhead Stirrers and Hei-Connect and Hei-End Magnetic Stirrers. The software allows to operate the device with all its functions entirely from a PC.

Version	Alteration date
2.0.0.4	12/2019

#### Variants

The software is available in different configurations. Certain features and functions apply to specific configurations only. The varieties are documented in this manual.

### > System requirements

System requirements	Minimal	Recommended
Operating system	Windows 7 (or higher) 32 Bit	Windows 7 (or higher) 64 Bit
Working memory RAM	2 GB	4 GB
Processor	2,5 GHz - Dual-Core-processor	2,5 GHz - Quad-Core-processor
Interface	1 x USB oder 1 x RS232	depending on amount of stirrers
Image	6.500 colors	16 million colors
Screen resolution	1400 x 1050 pixels	1680 x 1050 pixels

The software needs approx. 60 MB free storage space. If you would like to control more than one stirrer, your PC must meet the recommended system requirements.

### > About this manual

This operating manual is a component of the device described here. You will also find the manual on the USB stick in the directory: „program files\Heidolph\Hei-Control“.

- Please read this manual carefully and obey all safety and warning notices.
- Ensure that every operator reads this manual.
- Ensure that this manual is accessible for every operator.
- Pass on the operating manual to the subsequent owner.

For the current version of this operating manual in pdf format, please go to:

<http://www.heidolph.com/support/hei-control>

## ➤ Terms and signs used

In this manual you will find standardized terms and signs to warn you of possible dangers or give important hints. Please take special notice to these signs to avoid personal injury or damage to the appliance.

The following terms and type of signs (basic drawing elements) are used:

Symbol	Additional term / Description
<b>Warning signs</b> 	The yellow triangle indicates hazardous situations. It is used in combination with the following terms: <b>DANGER:</b> Indicates a hazardous situation which, if not avoided, <b>will</b> result in irreversible injury or death. <b>WARNING:</b> Indicates a hazardous situation which, if not avoided, could result in serious injury or material damage. <b>CAUTION:</b> Indicates a hazardous situation which, if not avoided, may result in a minor injury or material damage.
<b>Prohibitory signs</b> 	Not to do: The red circle indicates a situation that should be avoided under all circumstances to avoid injury or damage.
<b>Mandatory signs</b> 	To do: The blue circle indicates important information. Please obey to avoid property damage.

### Other signs used

Symbol	Description
→	Handling instruction, action required
✓	Result of action
▪ ▪ ▪	List of information
a. b. c.	List of variants

## ➤ **General safety instructions**

→ Please ensure the operating manual is available at all times.

### **The software may only be used under the following circumstances:**

- Only operate the software, if the herewith operated device is in full working order.
- Ensure all operators of the software possess the necessary safety and risk awareness for operating the device.
- If there is something you do not understand, or certain information is missing, ask your manager or contact the manufacturer.
- Do not operate the software without authorization.
- Only operate the software to use the device according to its intended use.

## **Intended use**

The software is intended for use by personnel only, trained and authorized to operate the stirrers Hei-TORQUE Precision, Hei-Connect and Hei-End.

The software is suitable for the following use:	<ul style="list-style-type: none"> <li>▪ Software based operating of the stirrers Hei-TORQUE Precision, Hei-Connect and Hei-End</li> <li>▪ Control up to 4 devices from one PC</li> </ul>
The operation of the software is suitable:	on any PC meeting the system requirements (see section "System requirements")

### **Improper use**

Any use which deviates from the software and device's intended use is considered to be improper. The manufacturer does not accept liability for any damages resulting from non-permitted uses. The risk is carried by the operator alone.

## ➤ **Operating company's obligations**

### **Changes to the software**

- No unauthorized changes may be made to the device.
- Unauthorized changes shall void the EC Declaration of Conformity and the device must not be operated any more.
- The manufacturer shall not be held liable for any damage, danger or injuries resulting from unauthorized changes or from operating the device other than as described in this manual.

### **Personnel safety**

→ Only qualified personnel who are authorized to work with Hei-TORQUE Precision, Hei-Connect and Hei-End may operate the software and the device.

### ➤ Safety during use

- Pay attention to all safety instructions specified in the manuals for the Overhead Stirrer Hei-TORQUE Precision, and the Magnetic Stirrers Hei-Connect and Hei-End while operating the unit through the software.
- Position the provided warning sign before you switch on the device.
- Do not start the device through the software when people are near to avoid any danger.
- Lock your PC to prevent unauthorized use when the device is linked to the PC.

### ➤ Disposal

- Dispose of the software carrier according to the valid national legal regulations.
- Dispose of the packaging material in accordance with the valid national legal regulations.

### ➤ Installation

Install the software before you connect the device with the PC.

The software is delivered on a USB stick.

- ➔ Insert the USB stick into the interface of your PC.

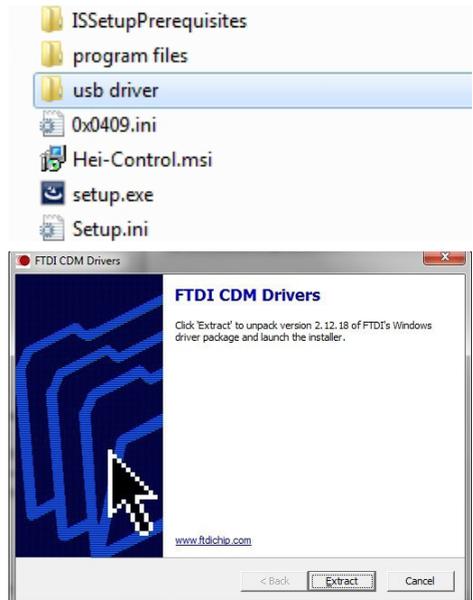
USB stick

### USB interface

To operate the software with a USB interface you must install an additional driver. The driver simulates a COM interface at the USB interface.


You can find the driver on the delivered USB stick.

- ➔ Insert the USB stick into the interface of your PC.
- ✓ Adjoining folders and files are shown on the removable medium.
- ➔ Open the folder „usb driver“.
- ➔ Double click „CDM v2.12.18 WHQL Certified“.
- ➔ Follow the instructions on the screen.
- ➔ Press „next“ to accept the storage location“.
- ✓ The driver is stored in the folder for Windows driver.
- ➔ Complete the installation with „Finish“.

### Software Hei-Control

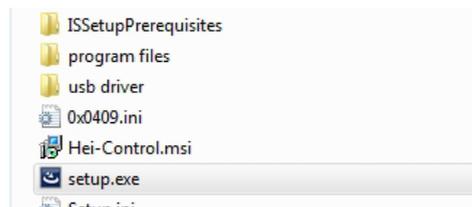- ➔ Insert the USB stick into the interface of your PC.
- ✓ Adjoining folders and files are shown on the removable medium.
- ➔ Double click „setup.exe“.


## Installation


If double click on „setup.exe“ does not work your system configuration requires a different method.

Please click „setup.exe“ with the right mouse button and execute the file as administrator.

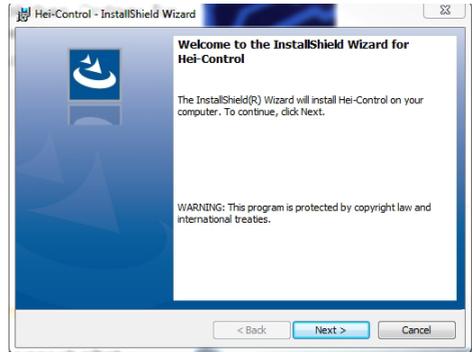- ✓ The start screen for installation opens.
- Click „Next“.
- Enter your user name via “user name“.
- Confirm with „Next“.

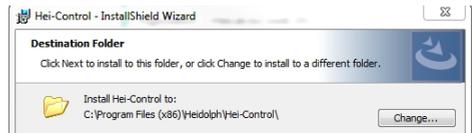- If applicable change the suggested storage location for the software Hei-Control via „Change“.
- Follow the instructions on the screen.
- Complete the Installation with „Finish“.

- ✓ The icon for executing the software is stored on your desktop.
- ✓ The exe-file is stored on the stated storage location.

## Manual Hei-Control

After completed installation, the manual “BAL Hei-Control EN-DE” is stored on the same storage location in which the exe-file is stored (see above).

You can also view the manual via the icon “Help” in the software menu.

# Installation adjustments

You can modify or repair the program installation through the program "setup.exe".

- Insert the USB stick into the interface of your PC.
- Double click "setup.exe".
- ✓ The program recognizes a pre-installed program-version and shows adjoining information.
- Click "Next".

- You receive adjoining options.

### Modify:

No function is provided with the free-software version.

### Repair:

The repair program replaces missing, accidentally deleted or shifted files. The program can't repair damaged files.

### Remove:

The remove program uninstalls the software Hei-Control.

- Choose the dedicated option.
- Follow the instructions on the screen.
- Complete the program with "Finish".

# Updates

If you receive an update for the current software version, the previous version must be deinstalled. You can uninstall the version through the appropriate windows program.

Saved profiles are still accessible.

## ➤ Interface

A serial interface is located on the back of the device.

### Interfaces Overhead Stirrers

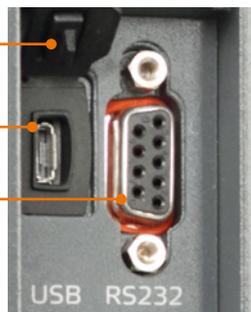
Depending on model their are two variants of interfaces:

- USB interface, Micro Type AB
- RS 232 interface, SUB D9

USB interface dust protection cap

USB interface

RS232 interface

### Interfaces Magnetic Stirrers

An RS 232 interface is located on the back of the device.

RS 232 interface

You can control up to 4 devices from one PC through several interfaces on the PC.

### Connect interfaces

- ➔ The device is switched off.
- ➔ Insert the plug of the delivered interface cable into the interface of the device.
- ➔ Insert the other end into the interface of your PC.

If your PC only disposes of USB interfaces, use a USB-RS 232 adaptor cable.

## Start software and link device

### Warning (Overhead Stirrer): Automatic start! Risk of injury! Risk of entanglement!

Rotating parts can injure the user and wind up loose ends of hair, clothing, jewelry and other items.

Keep loose hair, clothing and jewelry away from rotating parts that can harm the user. Position the provided warning sign before you activate remote-control. Every user must be aware of the potential risk of injury.

### Warning (Magnetic Stirrer): Risk of burning!

The hotplate can reach temperatures up to 300 °C.

If you work with temperatures above 50 °C wear safety gloves and appropriate protective clothing.

## Start and end program

### Start program

- Double click the icon "Hei-Control" .
- ✓ Adjoining start screen is shown.
- Select "Magnetic Stirrer" or "Overhead Stirrer".

Select your device

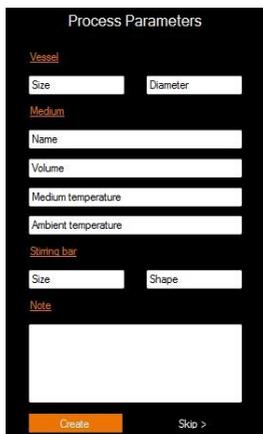
Magnetic stirrer

Overhead stirrer

- ✓ The "Process Parameters" window is displayed.
- Enter the parameters and click "Create" or "Skip" this step (see section "Process Parameters").
- The three-part menu of the software is displayed (see section "Menu structure")

### End program

- ✓ The stirring process is finished.
- Magnetic Stirrers:**
- ✓ The heating process is finished.
- Click the cross on the screen to end the program.

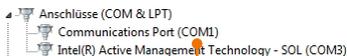
## Link device and PC

- Select "Ports" through the device manager of your PC.
- Check which COM ports are already occupied.

### Serial interface of PC

The serial interface is identified as "Communications Port" in the device manager.

That interface is always visible in the device manager, independent of the occupancy.

Physical serial interface

### USB interface of PC

The USB interface is only visible in the device manager when a USB cable is connected which simulates a serial interface.

Virtual serial interface

## Start-Up

- Plug in the interface cable into the device and PC.
- ✓ The COM interface identified as "serial" is shown in the device manager.
- Switch on the device.
- Start the software Hei-Control (see section "Start and end program").

- Choose the suitable COM interface through "Port" in the „Process setting“ area. If necessary, update the port list with „R“.

MR Hei-Connect

Port

- ✓ The connected device is listed. You can enter a dedicated name for the device.
- ✓ The display of the device blinks for a few seconds.
- Start the stirring process.

### For Magnetic Stirrers only:

- Start the heating process.
- ✓ Communication link between PC and device is completed.
- ✓ The device is switched to remote-operation.
- ✓ All commands are exclusively executed through the PC.

### Overhead Stirrer

- ✓ The slider switch of the stirrer is still active to end the stirring process.

### Magnetic Stirrer Hei-Connect

- ✓ The main functions can still be executed.
- ✓ You can always switch off the unit through the ON/OFF button.

If the software freezes while you link a device, check the driver installation of the USB interface and use the recommended USB-serial adaptor.

## Cut link between device and PC

- Stop the stirring process through the touch screen (Overhead Stirrers only) or using the ON/OFF button. The Magnetic Stirrer Hei-End must be unplugged.
- The communication link is cut off and the stirrer is manually operable.

If communication is disrupted, e.g. through disconnection of the interface cable, the device operates according the last settings for another 60 seconds and then comes to a complete stop. (Not valid for the magnetic Stirrer Hei-End.)

### ➤ Process parameters

After you started the program and selected the “Magnetic Stirrer”, the “Process Parameters” window is displayed (see section “Start Program”).

→ Enter the process parameters or skip this step with “Skip”.

#### Enter process parameters

→ Enter:

- size and diameter of the vessel
- name, volume and temperature of the medium
- ambient temperature
- type, size and material of the impeller
- a note

→ Click „Create”.

- ✓ The three-part software-menu is displayed (see section „Menu structure”).
- ✓ When you save process data (see section “Save and restore process data”), the process parameters are also saved in the .csv file.

**Process Parameters**

Vessel

Size Diameter

Medium

Name

Volume

Medium temperature

Ambient temperature

Impeller

Type

Size Material

Note

Create Skip >

#### View and correct process parameters

You can view and correct process parameters any time.

- Click the “Show process parameters” symbol  in the „Process setting” area.
- ✓ The “Process parameters” window is displayed.
- Adapt the parameters and click “Save”.

## ➤ Menu structure

The PC-surface enables you to view three different menus simultaneously, which are only available at the stirrer only via different sub-menus.

Via the enlargement icon  and downsizing icon  you can foreground one of the three menus as well as place the menus back.

### Main menu: Process setting:

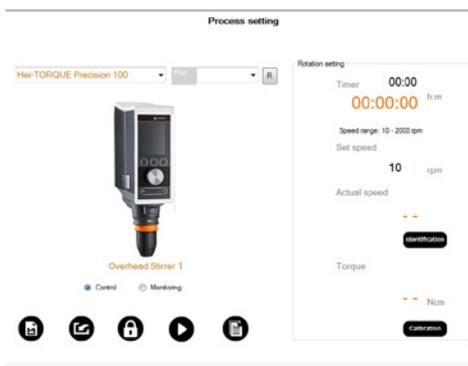
(see manual Hei-TORQUE, chapter "Device description, Hei-TORQUE Precision Display")

Here you choose the COM interface and the option for communication between your PC and the device.

With this menu you can enter stirring parameters and a stirring time.

The main menu is a reflection of the main menu on the device except "Menu" which is replaced by "Start".

- ➔ Use the cursor to move to entry fields for nominal values. Fields for actual values (orange characters) cannot be modified.
- ➔ Use the keyboard to type in numerical values.

Additionally, you can save and import process data, lock the software and view process parameters.


### Favorite menu: Programmable Control

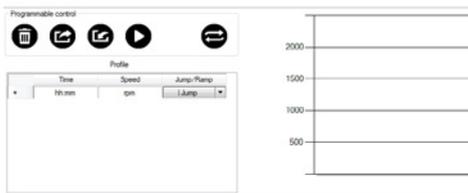(see manual Hei-TORQUE "Start-Up, Edit Profile")

Through the menu "Programmable Control" you can define profiles for a stirring process.

Move from field to field via cursor or keyboard.

- ➔ Use the cursor to click in the first field of the first row.
- ➔ Enter numerical values for "Time" and "Speed" through the keyboard.
- ➔ Choose the function "I Jump" or "/ Ramp".
- ➔ Confirm each row with Enter or Tab key.

# Overhead Stirrer Operation

## Graph menu

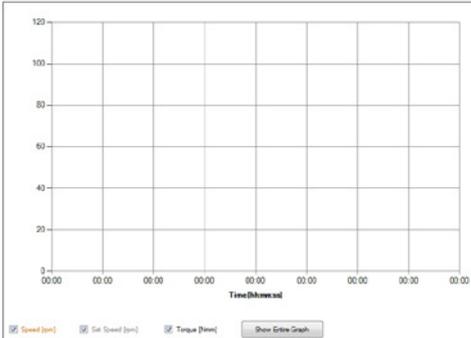
(see manual Hei-TORQUE, chapter "Operation, Display progression graph")

The process of a stirring operation is diagrammed. By default the last 5 minutes are shown.


Three factors can be displayed at the same time as well as separately:

- Actual speed
- Set speed
- Torque


→ Choose an option with the cursor.

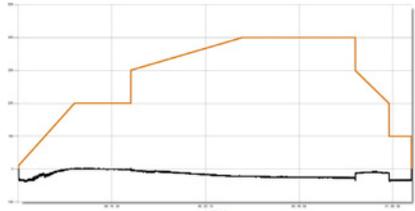
→ Choose "Show Entire Graph" to show the total process at this juncture.

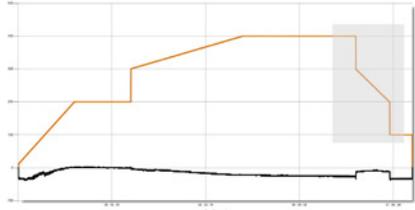→ Select the area you would like to view in detail.


✓ The area is enlarged.

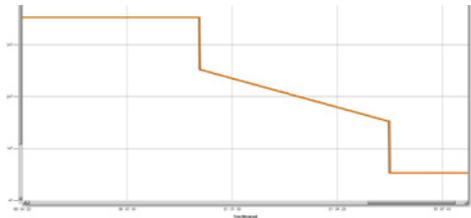→ Click on these symbols to undo zooming.

Undo vertical zoom

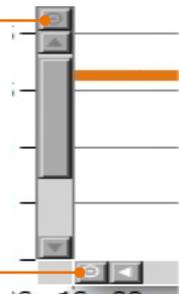Undo horizontal zoom

## Register and remove several Overhead Stirrers

You can register and control up to four devices from one PC.

### Register further stirrers

→ Click "New Device" and select "Overhead Stirrer".

- ✓ Another field for the next stirrer is integrated.
- To control the new integrated stirrer, choose the appropriate COM port (see section "Link device and PC").

### Remove stirrer

- ✓ The removable stirrer is not connected with the PC.
- Click the red cross next to the stirrer you want remove.
- ✓ The appropriate control panel disappears.

Closing the software removes all registered stirrers.

## Operate or monitor stirrer

In the "Process setting" area, you can choose from two options of communication.

### Operate stirrer

- Click "Control".
- ✓ The stirrer receives all operation commands through the software.

### Monitor stirrer

- Mark „Monitoring“.
- ✓ The stirrer can only be operated via device.
- ✓ The software reads the actual process.
- ✓ The process is displayed at your PC.

## Symbols used

### Area „Programmable control“

Symbol	Name	Function
	Clear	Reset profile
	Save	Save profile
	Import	Import saved profile
	Start	Start profile
	Pause	Interrupt stirring process
	Repeat	Repeat stirring process

### Area „Process setting“

	Save process	Save process data
	Restore process	Import process data
	Lock	Lock software
	Unlock	Unlock software
	Start	Start stirring process
	Stop	Stop stirring process
	Show process parameters	View and correct process parameters

- Black highlighted symbols (●) point to functions to choose of.
- Grayed out symbols (◐) point to functions which are not available and/or are currently activated.

## ➤ Control the Overhead Stirrer

### Set stirring parameters

- ✓ Stirrer and PC are connected with the interface cable.
- ✓ Stirrer and PC are switched on.
- Start the software Hei-Control.
- Set the favored stirrer and interface in the "Process setting" area.

#### Set range (Hei-TORQUE Precision 400 only)

- Go to the "Rotation setting" section of the "Process setting" area and click "Range.." to change gear.

#### Set speed

- Click in the box below "Set speed".
- Enter the dedicated speed through the keyboard.
- Confirm with return.

Set speed

100 rpm

- ✓ The set speed is shown below the box "Set speed".

Set speed  
120 rpm

#### Set Time (if required)

- Click in the box next to "Timer".
- Enter the dedicated time for the stirring process through the keyboard and confirm with return.

Timer 00:00  
00:00:00 h:m

## Start and Stop Stirring

**Warning:**  
**Automatic start!**  
**Risk of injury! Risk of entanglement!**

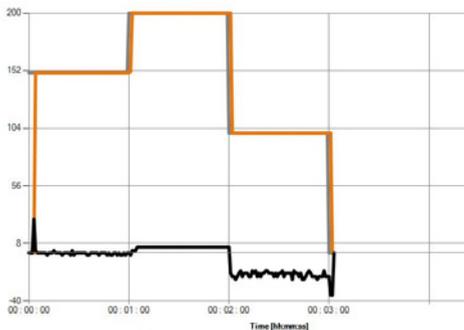
Rotating parts can injure the user and wind up loose ends of hair, clothing, jewelry and other items.

Keep loose hair, clothing and jewelry away from rotating parts that can harm the user.

Position the provided warning sign before you activate remote-control. Every user must be aware of the potential risk of injury.

### Start stirring

- Click the „Start“ symbol
- ✓ The stirring starts with the set speed.
- ✓ In the graph menu the stirring process is shown in a graph.
- ✓ The following factors are shown at the same time as well as separately:
  - Grey: Set speed
  - Orange: Actual speed
  - Black: Torque

### Stop rotation

- Click the „Stop“ symbol
- ✓ The stirring stops immediately.

You can also always stop the stirring on the device directly. There are two options:

- Stop the stirring through the slide bar.
- Switch off the device through the ON/OFF switch.

Both options stop the stirring process immediately.

## Profile

### Edit profile

- Maximum number of process steps: unlimited
- Maximum process time: 99 hours and 59 minutes

(See manual Hei-TORQUE, chapter "Edit profile")

- Enter the time of the first process step through the keyboard relative to zero.
- Enter the dedicated torque through "Speed".
- Choose from option Jump or Ramp through "Jump/Ramp".
- Complete each process step with return or tab.
- ✓ Every step of your input is shown graphically next to "Profile" on the right side.

Profile		
Time	Speed	Jump/Ramp
00:10	600	/ Ramp
00:20	600	/ Ramp
00:30	1200	/ Ramp
00:40	1200	Jump
00:50	1500	/ Ramp

### Save profile, recall and reset

#### Save profile

- Click the Save symbol
- Enter a name.
- ✓ The profile is saved under "My documents\Hei-Control".

#### Recall profile

- Click the Import symbol
- ✓ Previous saved profiles are shown.
- Double click onto the dedicated profile.
- ✓ The profile is imported with all edited standards and the graph diagram.

#### Reset Profile

- ✓ You have recalled a profile.
- Click the Clear symbol and confirm with "yes".
- ✓ Profile and graph are reset.
- ✓ You can edit and save a new profile.

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Save a new profile by using a previous profile as a template. Save the profile after editing under a new name.

# Overhead Stirrer Operation

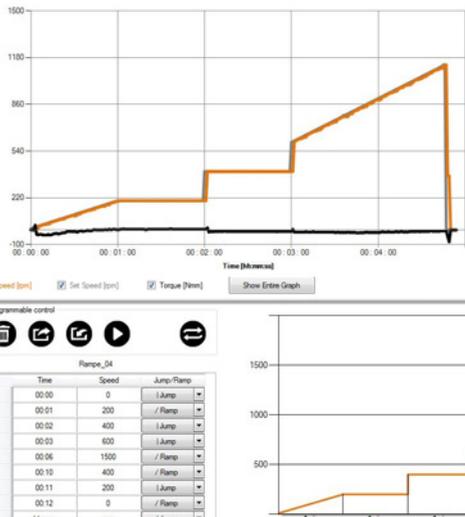
## Delete profile

- Open the register on your PC in which you have stored the profiles.
- Delete the profile not longer required.

## Start profile, interrupt, repeat, stop

### Start profile

- ✓ You have imported a profile.
- Click the Start symbol
- ✓ The stirring process starts.
- ✓ The process is shown, depending on the option, in the graph on the top left side.

### Interrupt profile

- Click the Pause symbol
- ✓ The process is halted.
- Click the Start symbol
- ✓ The stirring process continues from the halted position.

If the last process step of a profile does not end with speed „0“, the stirring process continues with the last values even it the time you set is over.

### Repeat profile

You can set the started profile into continuous operation.

- Click the Repeat symbol
- ✓ The color of the symbol switches from black to grey.
- ✓ The profile repeats until you stop the continuous operation.

To repeat a profile, the last process step in your profile must complete with torque "0".

The repetition can only start at standstill of the stirrer.

### Stop profile

- Click the Stop symbol
- ✓ The process stops.

### Calibrate torque

(See manual Hei-TORQUE, section "Calibrate torque")

- ✓ You have currently started stirring.
- In the "Rotation setting" section of the "Process setting" area, click on "Calibration".
- ✓ The torque is calibrated.
- ✓ In the graph diagram the torque goes to range negative and range positive.

Calibration

### Reset calibration

The calibration resets automatically when the stirring process stops.

### Read out Data

The function "Monitoring" allows you to operate the stirrer via device and at the same time shows you the graph of the stirring process on the PC.

- Mark "Monitoring" in the "Process setting" area.
- Start a stirring process on the device.
- ✓ The menu graph on the PC shows the stirring process.

Control  Monitoring

- Complete the stirring process.
- ✓ The graph of the stirring process stops.
- ✓ The communication to the PC is cut off.

The shown data will not be saved. As soon as a new stirring process is started or the program is shut down, the shown graph will be lost.

## Save and restore process data

### Save process data

- Click the „Save process" symbol .
- Enter a name.
- ✓ The data of all process steps are saved in .csv format under „My documents\Hei-Control".

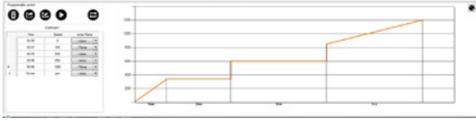
# Overhead Stirrer Operation

## Restore process data

- Click the "Restore process" symbol
- ✓ The register is opened in which you have stored previous process data.
- Double-click the dedicated process file.
- ✓ The process data is imported with all edited standards and the graph diagram.

Name	Änderungsdatum	Typ
Process1.csv	01.09.2016 14:34	Microsoft Office Excel-CSV
Process2.csv	24.08.2016 17:00	Microsoft Office Excel-CSV

When you open the .csv file in a text editor, the entered process parameters are displayed in addition to the table with the detailed process step data.

## Lock and unlock software

### Lock software

- Click the Lock symbol
- ✓ The dialog box „Password setting“ is displayed.
- ✓ Enter a password and confirm with „OK“.
- ✓ The input fields are locked and all symbols not to be used are not visible anymore.

### Unlock software

- Click the Unlock symbol
- In the dialog box „Password setting“, enter the password and confirm with „OK“.
- ✓ The input fields and all symbols are available again.

If you forget the password you entered, you must stop the device by closing the tab of the device.

## ➤ Process parameters

After you started the program and selected the "Magnetic Stirrer", the "Process Parameters" window is displayed (see section "Start Program").

→ Enter the process parameters or skip this step with "Skip".

### Enter process parameters

→ Enter:

- size and diameter of the vessel
- name, volume and temperature of the medium
- ambient temperature
- size and shape of the stirring bar
- a note

→ Click „Create“.

- ✓ The three-part software-menu is displayed (see section „Menu structure“).
- ✓ When you save process data (see section "Save and restore process data"), the process parameters are also saved in the .csv file.

**Process Parameters**

Vessel

Size  Diameter

Medium

Name

Volume

Medium temperature

Ambient temperature

Stirring bar

Size  Shape

Note

Create Skip >

### View and correct process parameters

You can view and correct process parameters any time.

- Click the "Show process parameters" symbol in the „Process setting“ area.
- ✓ The "Process parameters" window is displayed.
- Adapt the parameters and click "Save".

## ➤ Menu structure

The PC-surface enables you to view three different menus simultaneously, which are only available at the stirrer only via different sub-menus.

Via the enlargement icon  and downsizing icon  you can foreground one of the three menus as well as place the menus back.

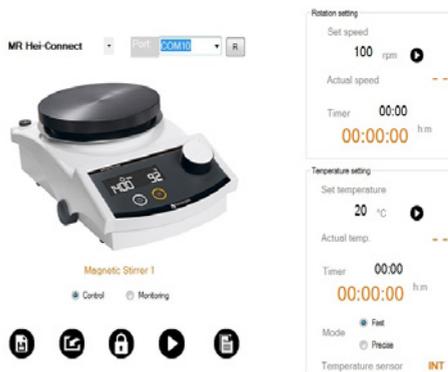
### Main menu: Process setting:

Here, you choose the COM interface and the option for communication between your PC and the device.

In this area you enter values for Speed, temperature and the duration of the stirring and heating process.

- Use the cursor to move to entry fields for nominal values. Fields for actual values (orange characters) cannot be modified.
- Use the keyboard to type in numerical values.

Additionally, you can save and import process data, lock the software and view process parameters.


### Favorite menu: Programmable control

Through the menu "Programmable control" you can define profiles for a stirring process.

Move from field to field via cursor or keyboard.

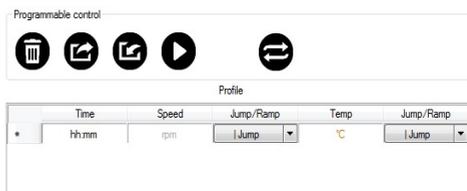- Use the cursor to click in the first field of the first row.
- Enter numerical values for "Time", "Speed" and "Temp" through the keyboard.
- Choose the function "I Jump" or "/" Ramp".
- Confirm each row with Enter or Tab key.

# Magnetic Stirrer Operation

## Graph menu

The process of stirring and heating operations is diagrammed. By default the last 5 minutes are shown.

Three factors can be displayed at the same time as well as separately:

- Actual speed
- Set speed
- Actual temperature
- Set temperature

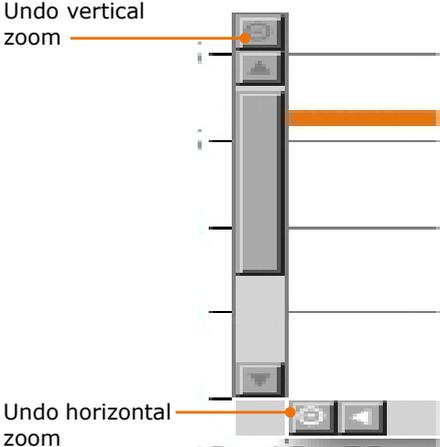
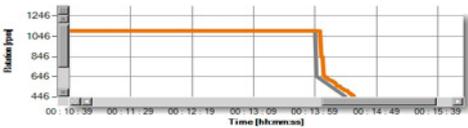
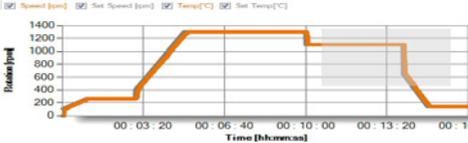
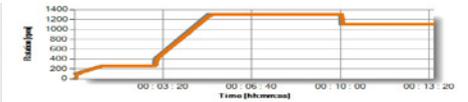
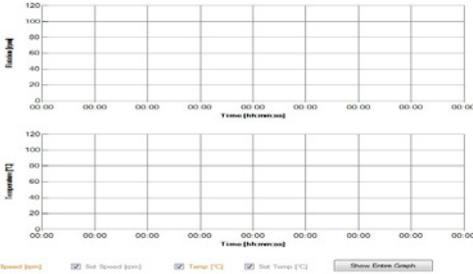
→ Choose an option with the cursor.

→ Choose "Show Entire Graph" to show the total process at this juncture.

→ Select the area you would like to view in detail.

✓ The area is enlarged.

→ Click on these symbols to undo zooming.

## Register and remove several Magnetic Stirrers

You can register and control up to four devices from one PC.

### Register further stirrers

→ Click "New Device" and select "Magnetic Stirrer" from the list.

- ✓ Another tab for the next stirrer is integrated.
- To control the new integrated stirrer, choose the appropriate COM port (see section "Link device and PC").

### Remove stirrer

- ✓ The Magnetic Stirrer to be removed is not connected with the PC.
- Click the red cross next to the stirrer you want to remove.
- ✓ The appropriate control panel disappears.

Closing the software removes all registered stirrers.

## Operate or monitor stirrers

In the "Process setting" area, you can choose from two options of communication.

### Operate the stirrer

- Click "Control".
- ✓ The stirrer receives all operation commands through the software.

### Monitor the stirrer

- Mark „Monitoring“.
- ✓ The stirrer can only be operated via device.
- ✓ The software reads the actual process.
- ✓ The process is displayed on the monitor of your PC.

## Symbols used

### Area „Programmable control“

Symbol	Name	Function
	Clear	Reset profile
	Save	Save profile
	Import	Import saved profile
	Start	Start profile
	Pause	Interrupt stirring and heating process
	Repeat	Repeat stirring and heating process

### Area „Process setting“

	Save process	Save process data
	Restore process	Import process data
	Lock	Lock software
	Unlock	Unlock software
	Start all	Start stirring and heating process
	Stop all	Stop stirring and heating process
	Show process parameters	View and correct process parameters

- Black highlighted symbols (🟩) point to functions to choose of.
- Grayed out symbols (🟨) point to functions which are not available and/or are currently activated.

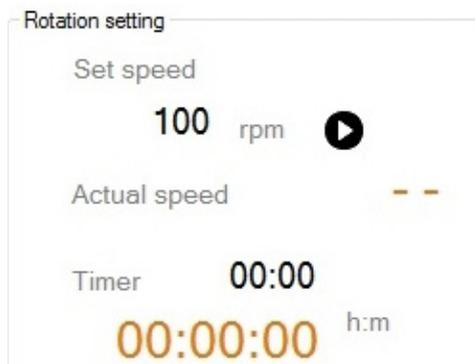
## ➤ Control the Magnetic Stirrer

- ✓ Stirrer and PC are connected with the interface cable.
- ✓ Stirrer and PC are switched on.
- Start the software Hei-Control.
- On the display of Hei-Connect, the blinking letters "PC" appear and on the display of Hei-End the word "Remote" is blinking.
- Set the favored stirrer and interface in the "Process setting" area.

## Set stirring parameters

You can set the parameters for stirring in the "Rotation setting" section of the "Process setting" area.

### Set speed

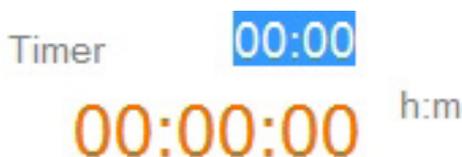
- Click the box below "Set speed".
- Enter the dedicated speed through the keyboard.
- Confirm with return.

- ✓ The set speed is shown below the box "Set speed".

### Set Time (if required)

- Click the box next to "Timer".
- Enter the dedicated time for the stirring process through the keyboard and confirm with return.


## Start and stop stirring

### Start stirring

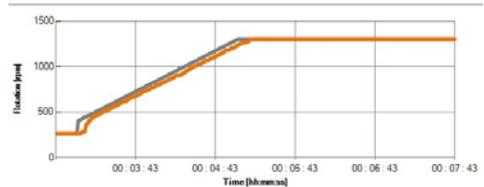- Click the „Start“ symbol  next to the value for the speed.
- ✓ The stirring starts with the set speed.
- ✓ In the graph menu, the stirring process is shown in a graph.

- ✓ The following factors are shown at the same time as well as separately:
  - Grey: Set speed
  - Orange: Actual speed

### Stop rotation

- Click the „Stop“ symbol  next to the value for the speed.
- ✓ The stirring stops immediately.

You can also always stop the stirring on the device directly. There are two options:

- Switch off the Magnetic Stirrer Hei-Connect through the ON/OFF switch.
- Unplug the Magnetic Stirrer Hei-End.

The stirring process stops immediately.

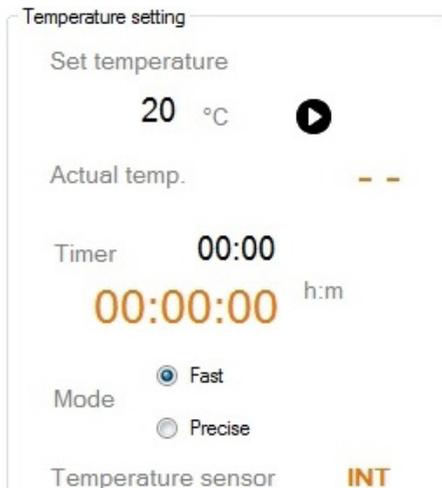
To start or stop the stirring and the heating process simultaneously, use the symbols „Start all“ and „Stop all“ in the lower part of the „Process setting“ area.

# Magnetic Stirrer Operation

## Set heating parameters

You can set the parameters for stirring in the "Temperature setting" section of the "Process setting" area.

### Set temperature

- Click in the box below "Set temperature".
  - Enter the dedicated temperature through the keyboard.
  - Confirm with return.
- ✓ The set temperature is shown below the box "Set speed".

Set temperature

20 °C

Set temperature

35 °C

### Set time (if required)

- Click in the box next to "Timer".
- Enter the dedicated time for the stirring process through the keyboard and confirm with return.

Timer 00:00  
00:00:00 h:m

As soon as the external temperature sensor Pt 1000 is connected, temperature will not be measured and controlled at the hotplate but directly at the sensor. The temperature displayed in the „Temperature setting” section is the temperature of the sensor and you can choose one of the modes: „Fast” or „Precise” (see manual Hei-Connect, Hei-End „Operation, External temperature sensor Pt 1000”).

## Start and stop heating

### **Warning:** **Risk of burning!**

The hotplate can reach temperatures up to 300 °C.

If you work with temperatures above 50 °C wear safety gloves and appropriate protective clothing.

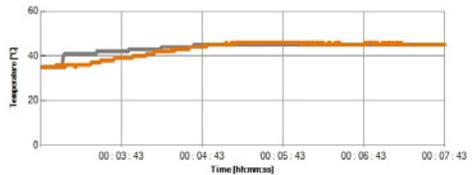
### **Warning:** **Risk of injury!**

Inflammable samples may ignite with high temperatures.

If you work with inflammable samples set temperature at least 25 °C lower than ignition point of sample.

### Start heating

- Click the „Start“ symbol  next to the value for the temperature.
- ✓ The heating starts with the set temperature.
- ✓ In the graph menu the heating process is shown in a graph.
- ✓ The following factors are shown at the same time as well as separately:
  - Grey: Set temperature
  - Orange: Actual temperature

### Stop heating

- Click the „Stop“ symbol  next to the value for the speed.
- ✓ The stirring stops immediately.

To start or stop the stirring and the heating process simultaneously, use the symbols „Start all“ and „Stop all“ in the lower part of the “Process setting” area.

## Profile

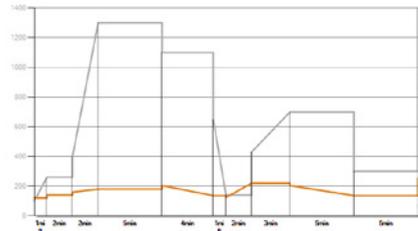
### Edit profile

- Maximum number of process steps: unlimited
- Maximum process time: 99 hours and 59 minutes

- Enter the time of the first process step through the keyboard relative to zero.
- Enter the dedicated "Speed".
- Enter the dedicated temperature in the "Temp" column.
- Choose from option Jump or Ramp through "Jump/Ramp".

test_1				
Time	Speed	Jump/Ramp	Temp	Jump/Ramp
00:00	100	I Jump	30	I Jump
00:01	250	/ Ramp	35	I Jump
00:03	400	I Jump	40	I Jump
00:05	1300	/ Ramp	45	/ Ramp
00:10	1100	I Jump	51	I Jump
00:14	650	I Jump	34	/ Ramp
00:15	140	/ Ramp	31	I Jump
00:17	430	I Jump	55	/ Ramp
00:20	700	/ Ramp	51	I Jump
00:25	300	I Jump	34	/ Ramp

- Complete each process step with return or tab.
- ✓ Every step of your input is shown graphically next to "Profile" on the right side.

If you enter a value for temperature and select the option „Jump“, the temperature does not reach the set value abruptly. The Magnetic Stirrer heats as quickly as possible with maximal power.

## Save profile, recall and reset

### Save profile

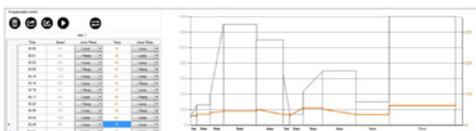
- Click the Save symbol
- Enter a name.
- ✓ The profile is saved under "My documents\Hei-Control".

### Recall profile

- Click the Import symbol
- ✓ Previously saved profiles are shown.
- Double click onto the dedicated profile.
- ✓ The profile is imported with all edited standards and the graph diagram.

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# Magnetic Stirrer Operation

## Reset Profile

- ✓ You have recalled a profile.
- Click the Clear symbol and confirm with "yes".
- ✓ Profile and graph are reset.
- ✓ You can edit and save a new profile.

Save a new profile by using a previous profile as a template. Save the profile after editing under a new name.

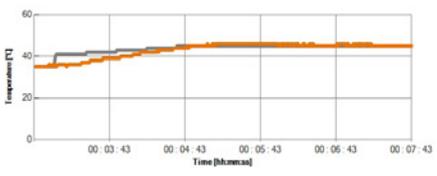
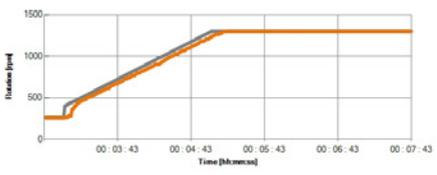
## Delete profile

- Open the register on your PC in which you have stored the profiles.
- Delete the profile not longer required.

## Start profile, interrupt, repeat, stop

### Start profile

- ✓ You have imported a profile.
- Click the Start symbol .
- ✓ The stirring and/or heating process starts.
- ✓ The process is shown, depending on the option, in the graph on the top left side.

### Interrupt profile

- Click the Pause symbol .
- ✓ The process is halted.
- Click the Start symbol .
- ✓ The stirring and heating process continue from the halted position.

Speed [rpm]  Set Speed [rpm]  Temp [°C]  Set Temp [°C]  Show Extra Graph

Programmable control

Time	Speed	Jump/Ramp	Temp	Jump/Ramp
00:00	150	/ Ramp	30	/ Jump
00:02	250	/ Ramp	35	/ Jump
00:03	400	/ Ramp	40	/ Jump
00:05	1200	/ Ramp	45	/ Ramp
00:10	1500	/ Ramp	51	/ Jump
00:14	600	/ Ramp	34	/ Ramp
00:15	150	/ Ramp	31	/ Jump

If the last process step of a profile does not end with speed and temperature „0“, the stirring process continues with the last values even it the time you set is over.

## Magnetic Stirrer Operation

### Repeat profile

You can set the started profile into continuous operation.

- Click the Repeat symbol .
- ✓ The color of the symbol switches from black to grey.
- ✓ The profile repeats until you stop the continuous operation.

To repeat a profile, the last process step in your profile must complete with torque "0".

The repetition can only start at standstill of the stirrer.

### Stop profile

- Click the Stop symbol .
- ✓ The process stops.

### Read out Data

The function "Monitoring" allows you to operate the stirrer via device and at the same time shows you the graph of the stirring process on the PC.

- Mark "Monitoring" in the "Process setting" area.
- Start a stirring or heating process on the device.
- ✓ The menu graph on the PC shows the stirring and heating process.
- Complete the stirring and heating processes.
- ✓ The graph of the stirring and heating processes stops.
- ✓ The communication to the PC is cut off

Control

Monitoring

The shown data will not be saved. As soon as a new stirring process is started or the program is shut down, the shown graph will be lost.

## Save and restore process data

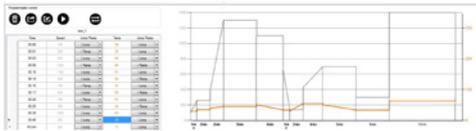
### Save process data

- Click the „Save process“ symbol
- Enter a name.
- ✓ The data of all process steps are saved in .csv format under „My documents\Hei-Control“.

### Restore process data

- Click the „Restore process“ symbol
- ✓ The register is opened in which you have stored previous process data.
- Double-click the dedicated process file.
- ✓ The process data is imported with all edited standards and the graph diagram.

Name	Änderungsdatum	Typ
Process1.csv	01.09.2016 14:34	Microsoft Office Excel-CSV
Process2.csv	24.08.2016 17:00	Microsoft Office Excel-CSV


When you open the .csv file in a text editor, the entered process parameters are displayed in addition to the table with the detailed process step data.

## Lock and unlock software
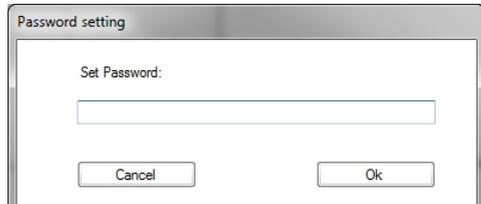
### Lock software

- Click the Lock symbol
- ✓ The dialog box „Password setting“ is displayed.
- ✓ Enter a password and confirm with „OK“.
- ✓ The input fields are locked and all symbols not to be used are not visible anymore.

### Unlock software

- Click the Unlock symbol
- In the dialog box „Password setting“, enter the password and confirm with „OK“.
- ✓ The input fields and all symbols are available again.

If you forget the password you entered, you must stop the device by closing the tab of the device.

## ➤ Malfunctions and clear up

General error messages	Possible reason	Troubleshooting
Do not change while in process!	Attempt to change parameters during operation	Stop operation, change parameters
The program has been modified or is already running!	exe file was renamed	Rename EXE file to original
	Program already running	Confirm with OK
Connection error!	<ul style="list-style-type: none"> <li>Interface cable not plugged in</li> <li>Wrong COM port assigned</li> </ul>	<ul style="list-style-type: none"> <li>Check plugs</li> <li>Change COM port (see „Stirrer parameters“)</li> </ul>
No data to save	No data was entered under "Profile"	Enter Data
Impossible to save profile	Save profile was cancelled	Save profile with "Save"
Impossible to restore profile: Wrong parameters in file	Profile contains invalid data	Enter a new Profile and save
Impossible to restore profile	Undefined error while importing profile	Enter Profile anew and save
Port xxx can not be opened!	COM port is used by other device	Change COM port

Hei-TORQUE Precision error messages	Possible reason	Troubleshooting
<ul style="list-style-type: none"> <li>Don't change while in process!</li> <li>Impossible to switch range!</li> </ul>	<p><b>Only Hei-TORQUE Precision 400:</b></p> <p>Attempt to change parameters during operation</p>	Stop operation, change parameters
The stirrer type is false! The connected stirrer is...	Wrong stirrer type entry	Enter correct stirrer in the area "process setting"
Error: Overload	<p>Motor or power adaptor overload:</p> <ul style="list-style-type: none"> <li>torque too high</li> <li>impeller blocked</li> </ul>	<p>Rotation stops automatically:</p> <ul style="list-style-type: none"> <li>remove cause for blocking,</li> <li>re-activate rotation</li> </ul>
Impossible to configure the device!	Stirrer could not be identified	Connected stirrer is possibly not a Heidolph unit

## Troubleshooting

<b>Hei-Connect, Hei-End error messages</b>	<b>Possible reason</b>	<b>Troubleshooting</b>
Hot plate sensor shorted!	Short circuit of hot plate sensor	Service matter; please contact your lab supplies distributor.
Hot plate sensor break!	Fracture of hot plate sensor	Service matter; please contact your lab supplies distributor.
Hot plate safety sensor shorted!	Short circuit of hot plate safety sensor	Service matter; please contact your lab supplies distributor.
Hot plate safety sensor break!	Fracture of hot plate safety sensor	Service matter; please contact your lab supplies distributor
Hot plate sensor difference >15°C!	Hot plate sensor; difference between sensors > 15°C	Service matter; please contact your lab supplies distributor.
Set safety $\Delta T$	Safety temperature $\Delta T$ exceeded; Indication of exothermic reaction: Controlling of temperature too slow	<ul style="list-style-type: none"> <li>Perform reset (power switch off/on)</li> <li>Increase value for safety temperature <math>\Delta T</math>, if applicable.</li> </ul> Service matter; please contact your lab supplies distributor.
External sensor shorted!	Short circuit of external temperature sensor	<ul style="list-style-type: none"> <li>Replace external temperature sensor</li> <li>Perform reset (power switch off/on)</li> </ul>
External sensor break!	Removal or fracture of external safety temperature sensor	<ul style="list-style-type: none"> <li>Insert or replace external temperature sensor</li> <li>Perform reset (power switch off/on)</li> </ul>
External safety sensor shorted!	Short circuit of external safety temperature sensor	<ul style="list-style-type: none"> <li>Replace external temperature sensor</li> <li>Perform reset (power switch off/on)</li> </ul>
External safety sensor break!	Removal or fracture of external safety temperature sensor	<ul style="list-style-type: none"> <li>Insert or replace external temperature sensor</li> <li>Perform reset (power switch off/on)</li> </ul>
External sensor difference >15°C!	External temperature sensor; difference between sensors > 15°C	<ul style="list-style-type: none"> <li>Replace external temperature sensor</li> <li>Perform reset (power switch off/on)</li> </ul>
Communication fault!	Internal communication error of device	Service matter; please contact your lab supplies distributor.
Triac error!	Defect of TriAC component	Service matter; please contact your lab supplies distributor.
Motor error!	Motor defect	Service matter; please contact your lab supplies distributor.
Kernel error!	Kernel defect (operating system)	Service matter; please contact your lab supplies distributor.

## Troubleshooting

The unit is stopped manually!	Heating or rotation switched off manually. (Controlling via PC is blocked.)	Restart device using PC software. (Blocking is switched off automatically.)
Hot plate sensor break or shorted (by hardware)!	Fracture or short circuit of hot plate sensors	Service matter; please contact your lab supplies distributor.
Hot plate safety $\Delta T$ exceeded!	Hot plate temperature has exceeded safety temperature $\Delta T$ ; indication of exothermic reaction: controlling of temperature too slow	Perform reset (power switch off/on) Service matter; please contact your lab supplies distributor.
External probe outside media!	External temperature sensor not inserted into sample	<ul style="list-style-type: none"> <li>▪ Insert the external temperature sensor into sample</li> <li>▪ And/or position container on the hot plate</li> <li>▪ Perform a reset (power switch off/on).</li> </ul>
External probe outside media (by start)!	At start of heating function, external temperature sensor not inserted into sample	<ul style="list-style-type: none"> <li>▪ Insert the external temperature sensor into the sample</li> <li>▪ And/or position container on the hot plate</li> <li>▪ Perform a reset (power switch off/on).</li> </ul>
External safety $\Delta T$ exceeded!	Safety temperature $\Delta T$ of the external temperature sensor exceeded; indication of exothermic reaction: controlling of temperature too slow:	<ul style="list-style-type: none"> <li>▪ Perform a reset (power switch off/on)</li> <li>▪ Service matter; please contact your lab supplies distributor</li> </ul>
Button heating!	Heating button defective	Service matter; please contact your lab supplies distributor.
Button motor!	Rotation button defective	Service matter; please contact your lab supplies distributor.
Button encoder!	Select knob defective	Service matter; please contact your lab supplies distributor.
Motor error!	Motor defective	Service matter; please contact your lab supplies distributor.
External sensor state changed during heating active!	Status external temperature sensor was changed during heating operation (while heating function was active)	Status external temperature sensor was changed during heating operation (while heating function was active)

If you experience a malfunction which can not be resolved, please contact your authorized Heidolph distributor immediately.

## ➤ Contact / Technical Service

### Questions / Repair work

If any aspect of installation or operation remains unanswered in the present manual, please get in touch with the following address or your local authorized Heidolph distributor.

#### Contact details:

---

#### **Heidolph Instruments Germany**

Heidolph Instruments GmbH & Co. KG  
Jürgen Heyder  
Walpersdorfer Str. 12  
D-91126 Schwabach / Germany

Phone: +49 - 9122 - 9920-13  
E-Mail: [juergen.heyder@heidolph.de](mailto:juergen.heyder@heidolph.de)

[www.heidolph.de](http://www.heidolph.de)

#### **Heidolph Instruments North America**

Phone: 1-866-650-9604  
E-mail: [service@heidolph.com](mailto:service@heidolph.com)  
[www.heidolphNA.com](http://www.heidolphNA.com)

#### **Heidolph Instruments United Kingdom**

Phone: 01799 - 5133-20  
E-mail: [service@radleys.co.uk](mailto:service@radleys.co.uk)  
[www.heidolph-instruments.co.uk](http://www.heidolph-instruments.co.uk)

#### **All other countries**

You will find contact details of your local Heidolph distributor at [www.heidolph.com](http://www.heidolph.com)

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This is a translation of the original operating manual. We reserve the right to make any technical modifications without prior notification. This documentation is not subject to revision; the latest version of this documentation can be found on our homepage. The copyright for all texts and images is held by Heidolph Instruments GmbH & Co. KG.
© Heidolph Instruments GmbH & Co. KG